

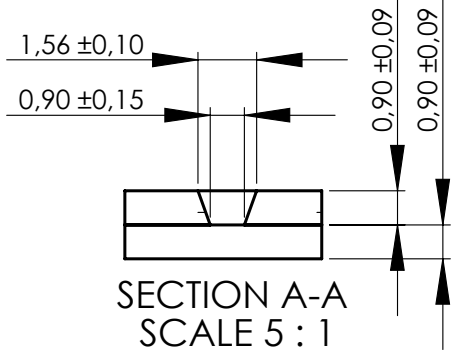
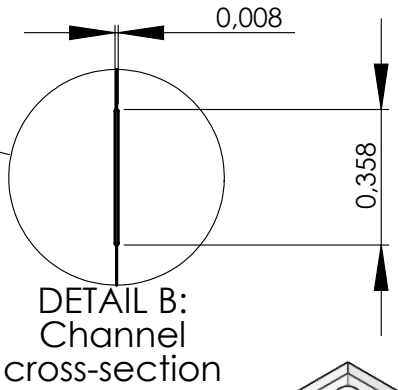
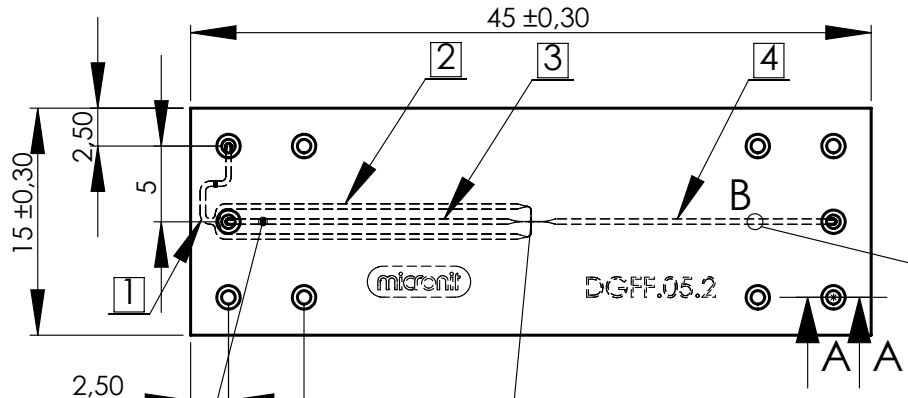



		Length [mm]	#	Volume [μ l]
1	Continuous inlet channel	7.75	1	0.023
2	Continuous channel	20.8	2	0.062 (0.124 tot.)
3	Dispersed channel	18.5	1	0.055
4	Outlet channel	18.5	1	0.055

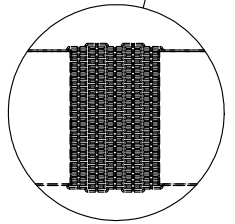

Nozzle dimensions:
0.005 x 0.0072

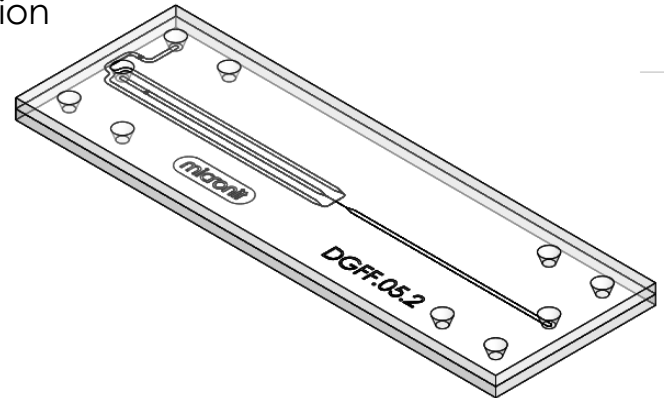

Filter structure at continuous inlet channel and dispersed channel

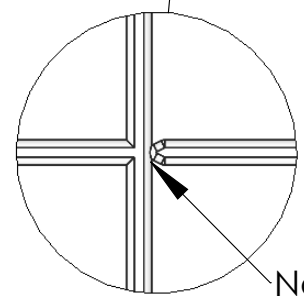

DETAIL C
SCALE 250 : 1

<p>THIRD ANGLE PROJECTION</p>	UNLESS OTHERWISE SPECIFIED: DIMENSIONS ARE IN MILLIMETERS			SURFACE FINISH: -	Sheet format: A4
	SURFACE TREATMENT: -			TOLERANCES: LINEAR: ±. ANGULAR: ±.	DEBUR AND BREAK SHARP EDGES
DRAWN	BvD	SIGNATURE	DATE 31-3-2017	TITLE: DGFF.05.2	
MOD					
CHK'D	WB		3-4-2017		
PROPRIETARY AND CONFIDENTIAL THE INFORMATION CONTAINED IN THIS DRAWING IS THE SOLE PROPERTY OF MICRONIT MICROFLUIDICS BV. ANY REPRODUCTION IN PART OR AS A WHOLE WITHOUT THE WRITTEN PERMISSION OF MICRONIT MICROFLUIDICS BV IS PROHIBITED.		Projectnr.	RD0961.002	DWG NO.	DGFF.05.2 R1
		MATERIAL:	Borofloat Glass	SCALE: 2:1	DO NOT SCALE DRAWING
		WEIGHT: 2.96			REVISION 1
					SHEET 1 OF 1