

micronit




Fluidic Connect Pro


quick start guide v1.4

micronit.com



0)

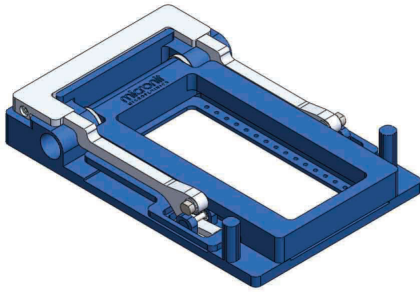


8x

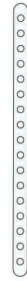

2x

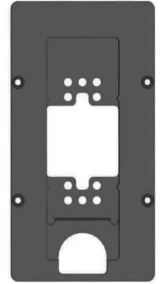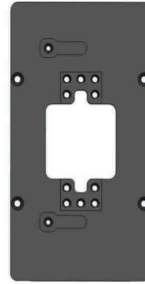

Part list

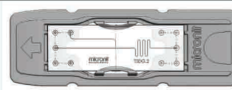

Example chip with connections
(sold separately)




3x




3x

1)

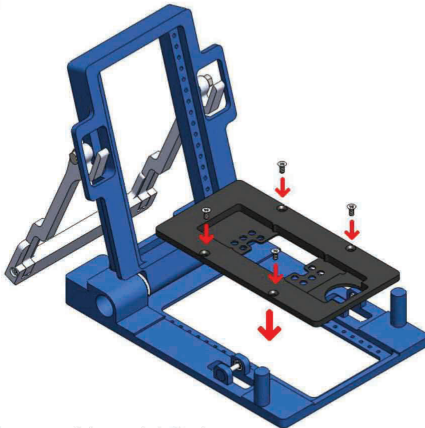


Place and (gently) fix bottom insert

2a) *optional for chips with thicknesses 0.9-1.4mm*

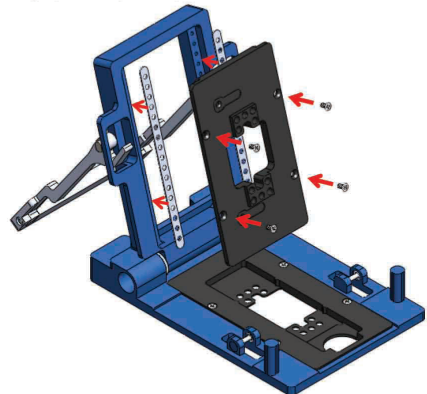


Place and fix spacers and top insert

2b)

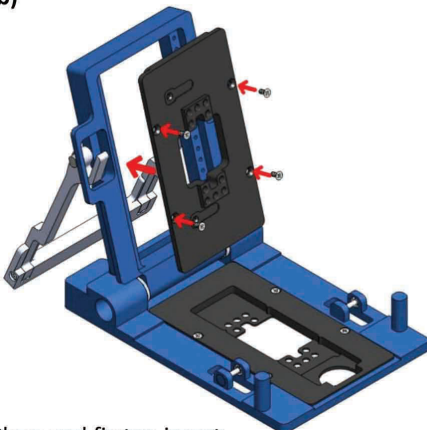


Place and fix top insert

3)

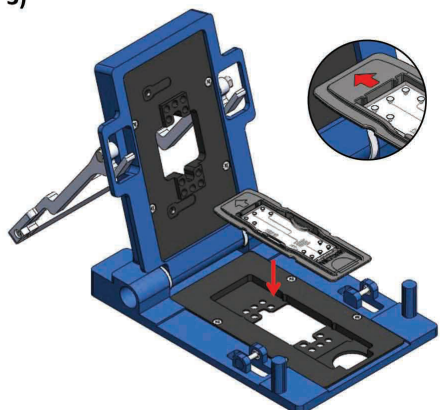


Position Fluidic Slide with microfluidic chip

4)

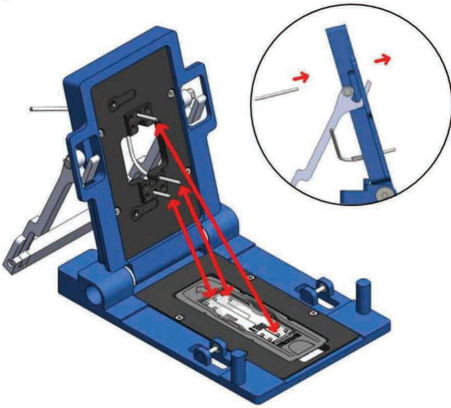


Slide the tubing in from the top side

5)

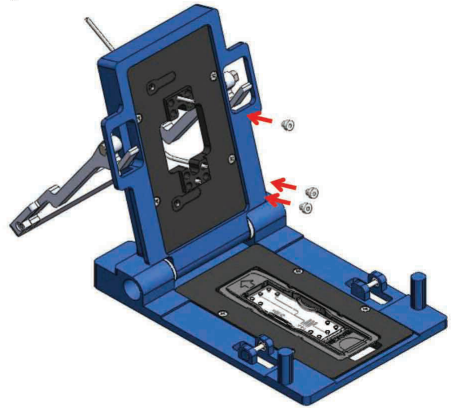


Assemble ferrules onto the tube ends

6)

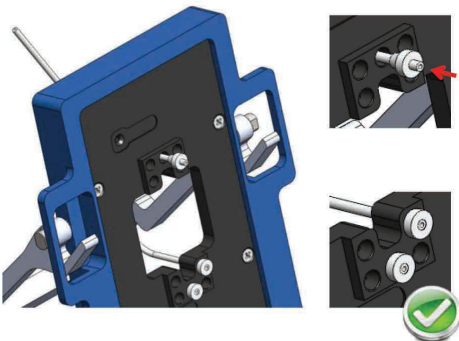


Press the tube with ferrule into the top insert

7)

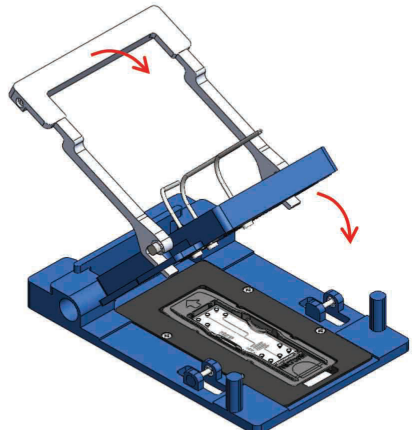


Close the holder using the handle

8)

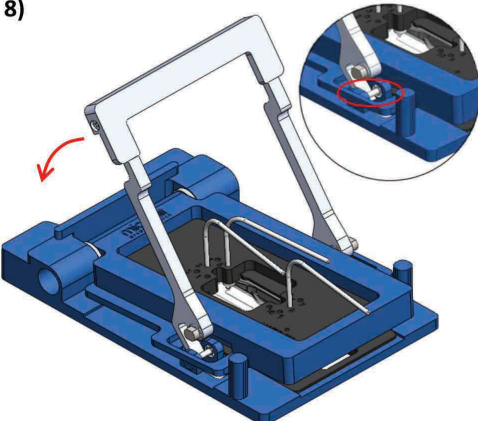

Fix the assembly using the gripper

9)

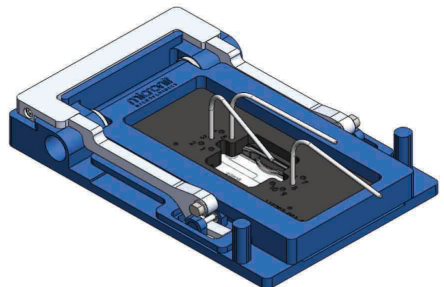

Your Fluidic Connect Pro is ready to use



Visit micronit.com for more information
about our products and services

