



# micronit

M I C R O F L U I D I C S


## Fluidic Connect Pro Resealable 4515 Quick Start Guide v2.0

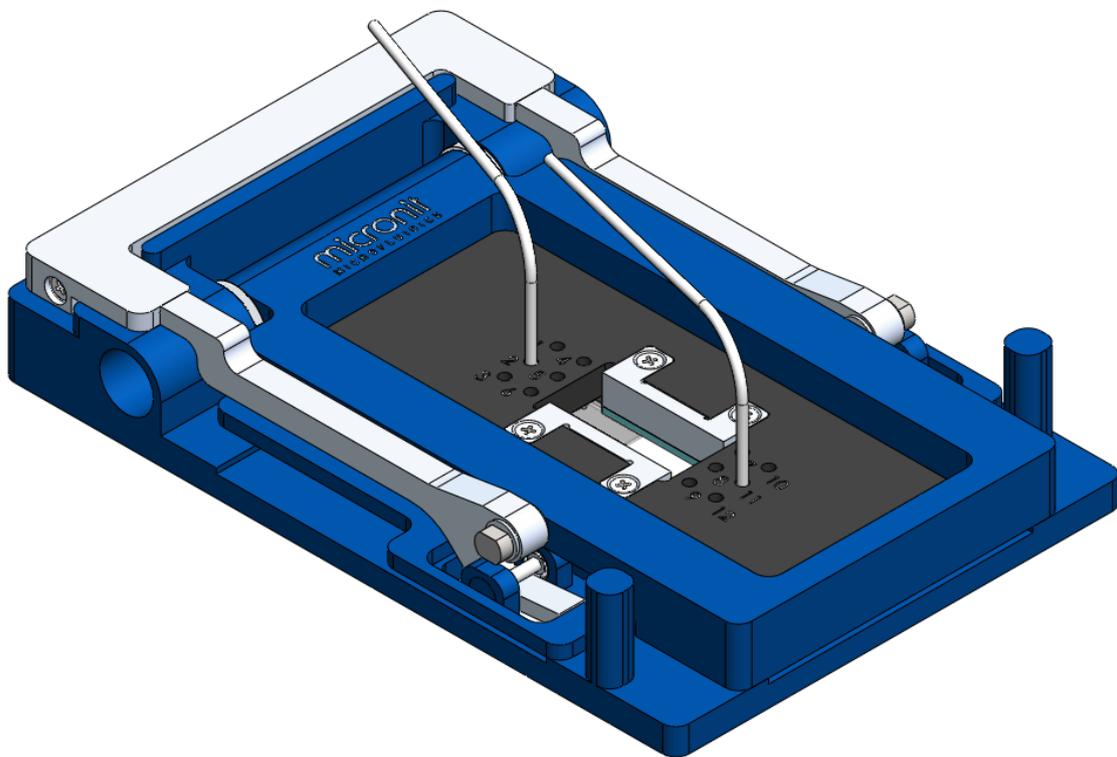



## Assembling the Fluidic Connect Pro – Resealable flowcell 4515

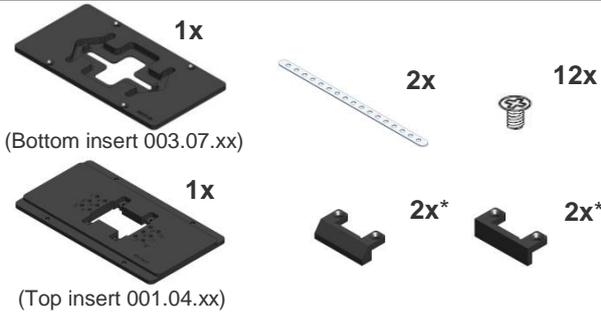

### For straight resealable Flowcell

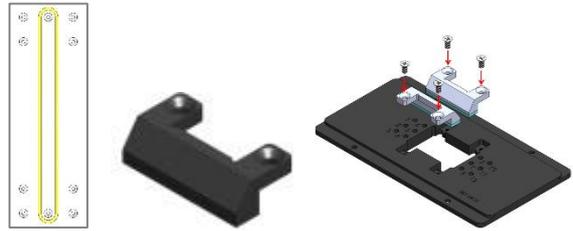


### For extended / wide resealable Flowcell

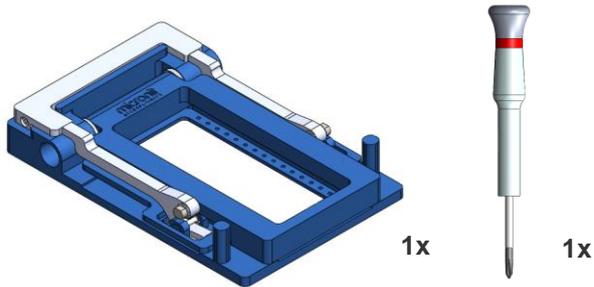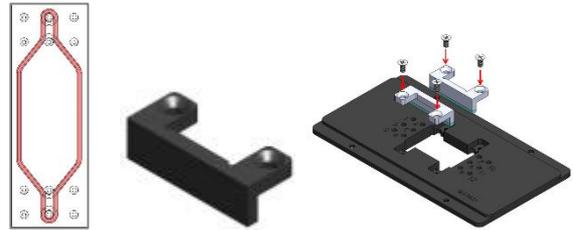


**1)** Be sure that the correct press inserts are mounted in the top insert.


\*Depending on Flowcell type (see step 1)

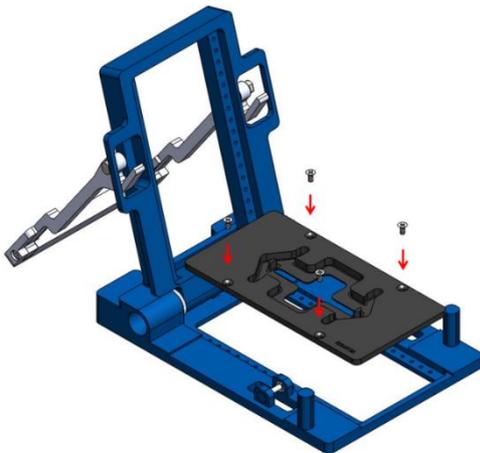


**2)** Place and fix bottom insert.

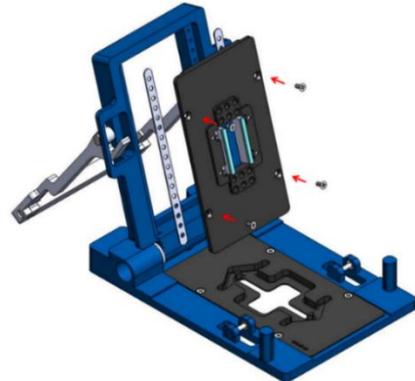

**3)** Place and fix top insert. Use the 0.5mm spacers between the top insert and Fluidic Connect Pro.






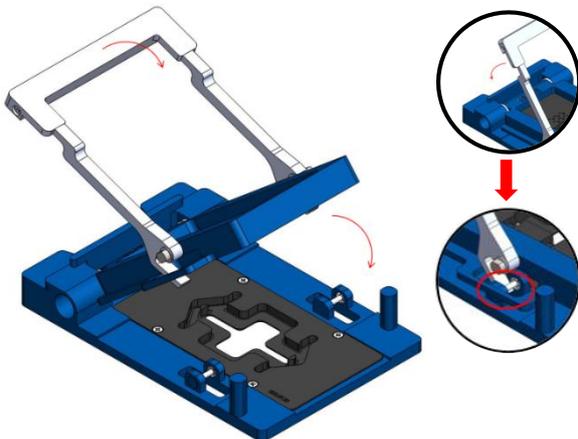


**4)** Close the holder using the handle and fix the assembly using the gripper.

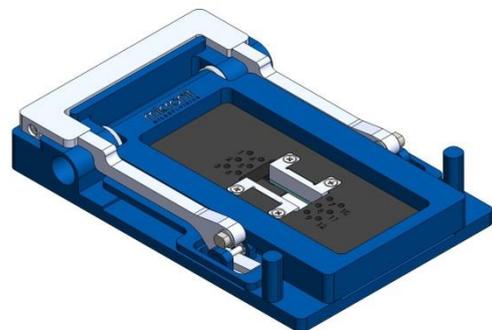

**5)** Your Fluidic Connect Pro is ready to use. (Assemble the fittings before making fluidic connections)

## Assembly of the fittings

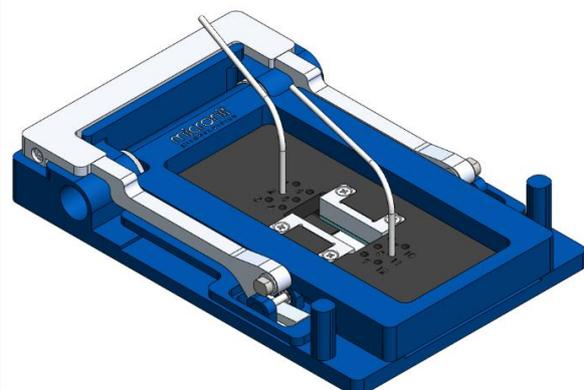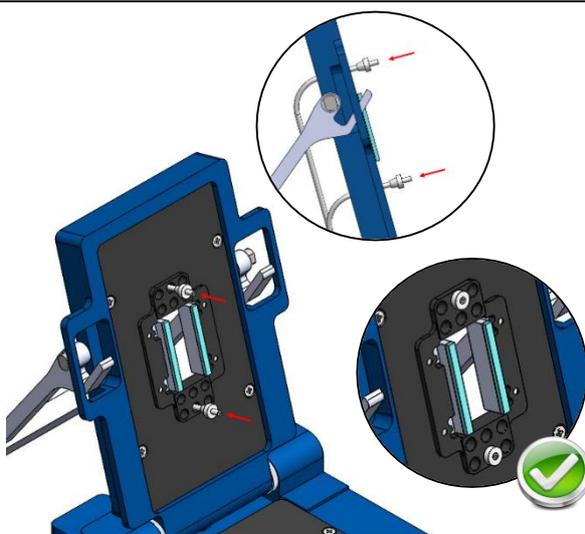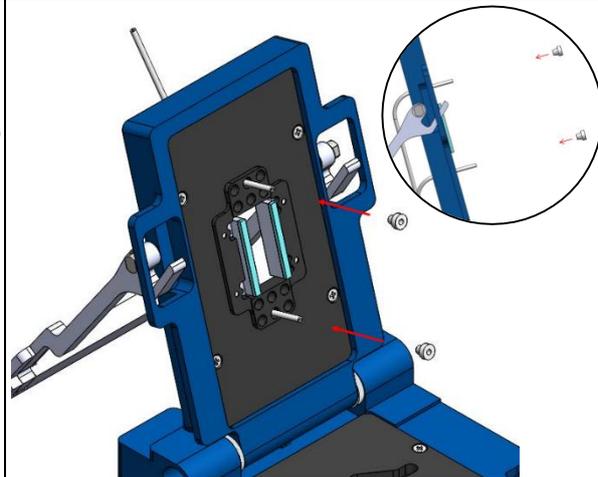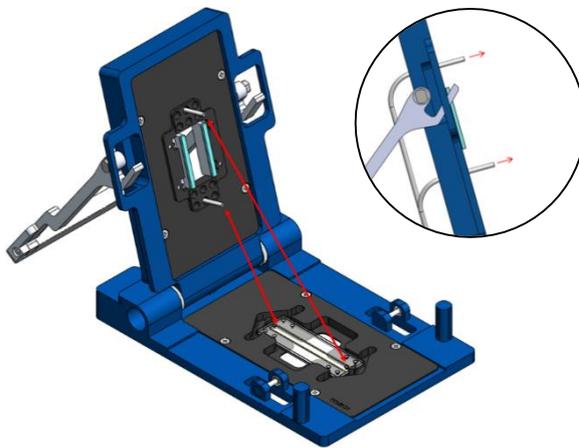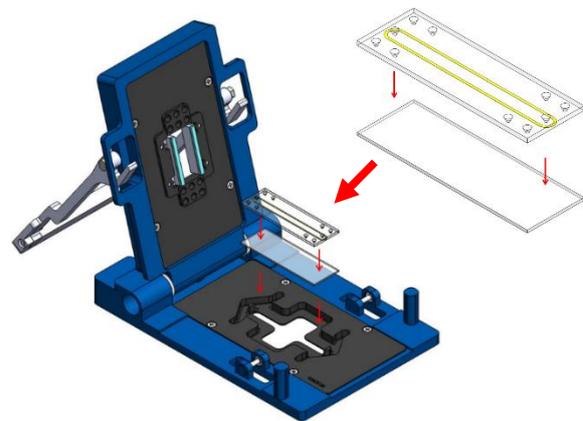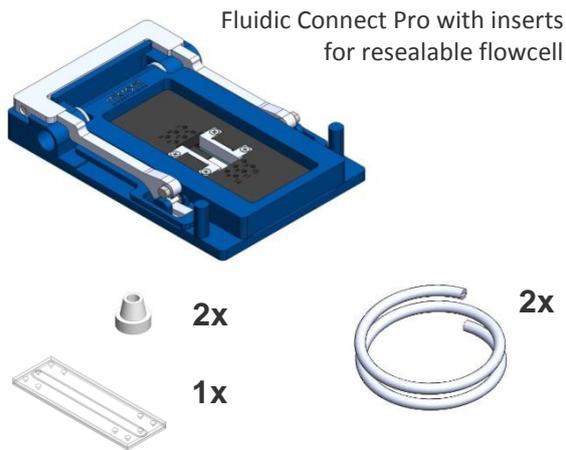

For more information go to: <http://store.micronit.com/chipholder/fluidic-connect-pro.html>

**micronit**  
MICROFLUIDICS

Address Colosseum 15 · 7521 PV Enschede · The Netherlands

Phone +31 53 850 6 850 Fax +31 83 850 6 851 E-mail [info@micronit.com](mailto:info@micronit.com) Internet [www.micronit.com](http://www.micronit.com)

